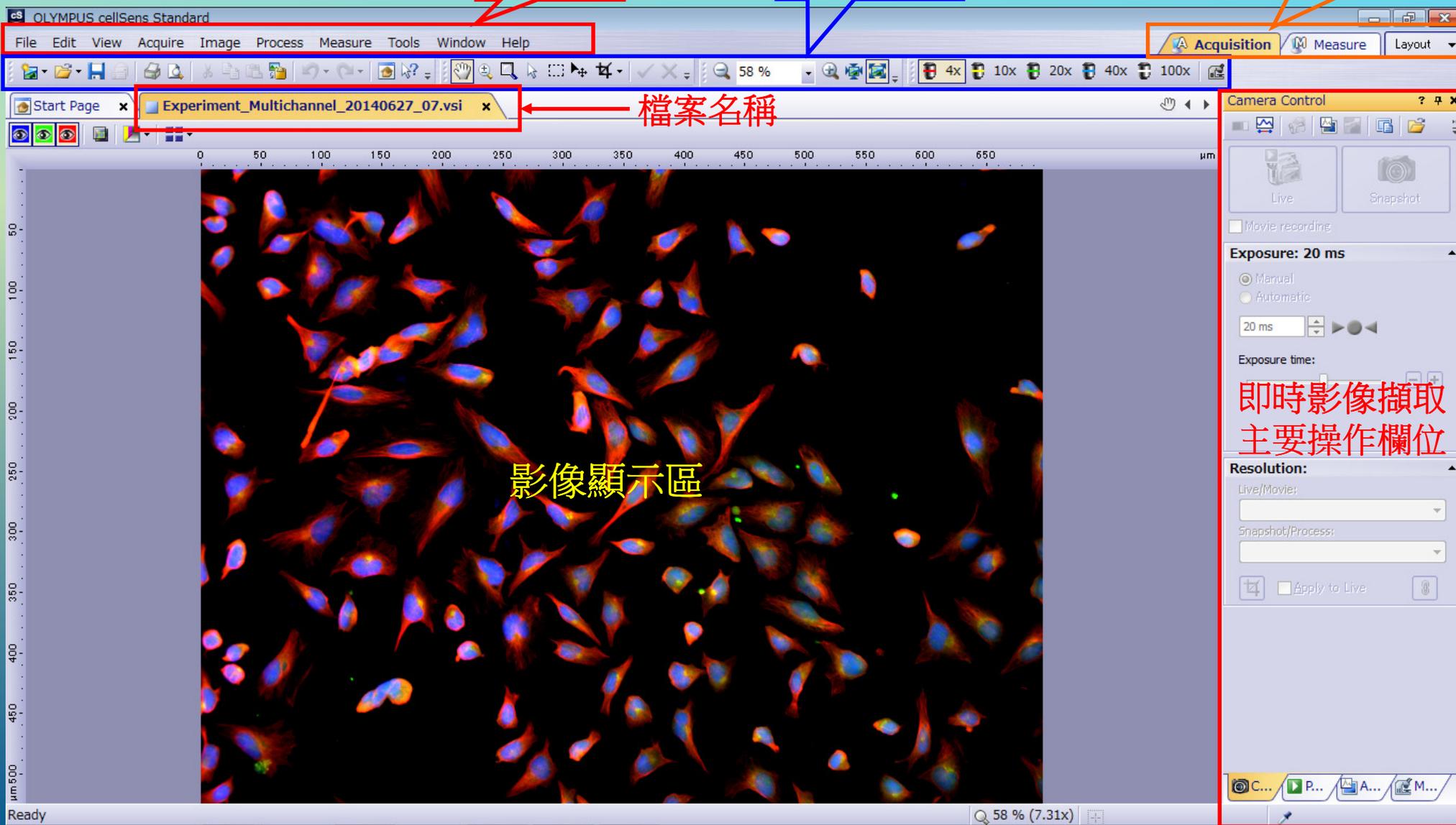
OLYMPUS cellSens
簡易操作

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① 開啟 cellSens

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
The screenshot shows the 'Camera Control' software interface with several settings and annotations:

- 彩色 黑白模式切換鍵 (部分CCD無此功能)**: Points to the color/black mode toggle button.
- 拍照擷取儲存設定**: Points to the snapshot/live recording buttons.
- Adjust Display**: Points to the display adjustment button.
- 白平衡 明視野使用**: Points to the white balance setting.
- 黑平衡 螢光使用**: Points to the black balance setting.
- 對焦輔助**: Points to the focus assist button.
- ⑥**: Points to the 'Movie recording' checkbox.
- ①**: Points to the 'Manual' exposure mode selection.
- ②**: Points to the 'SFL' (Spot Focus Lock) button.
- ④**: Points to the exposure time lock button.
- ⑤**: Points to the exposure range lock button.
- ⑦**: Points to the 'ISO 200' sensitivity setting.
- ③**: Points to the 'Exposure compensation' setting.
- ⑧**: Points to the 'Resolution: 1024 x 800' setting for Live/Movie.
- ⑨**: Points to the 'Resolution: 1024 x 800' setting for Snapshot/Process.
- 即時影像裁切**: Points to the crop button for live video.
- 動靜態畫素統一**: Points to the 'Apply to Live' checkbox.
- 曝光範圍置中**: Points to the exposure range lock button.
- SFL螢光專用模式(拍螢光時請打勾)某些CCD沒有**: Points to the SFL button.

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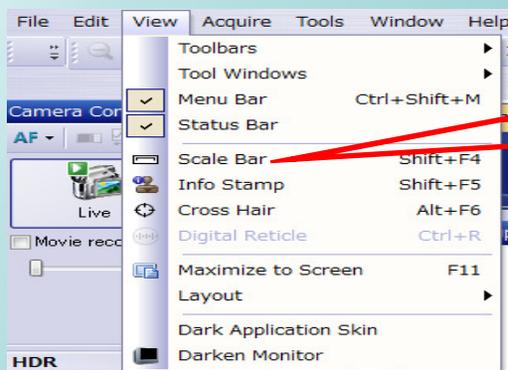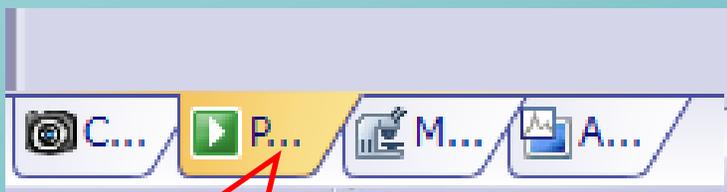
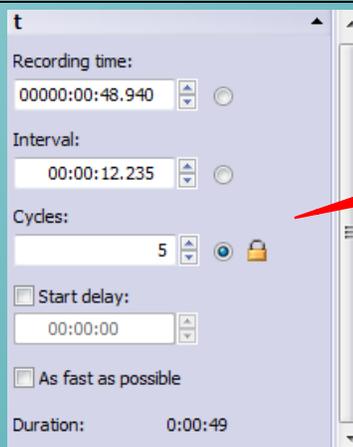
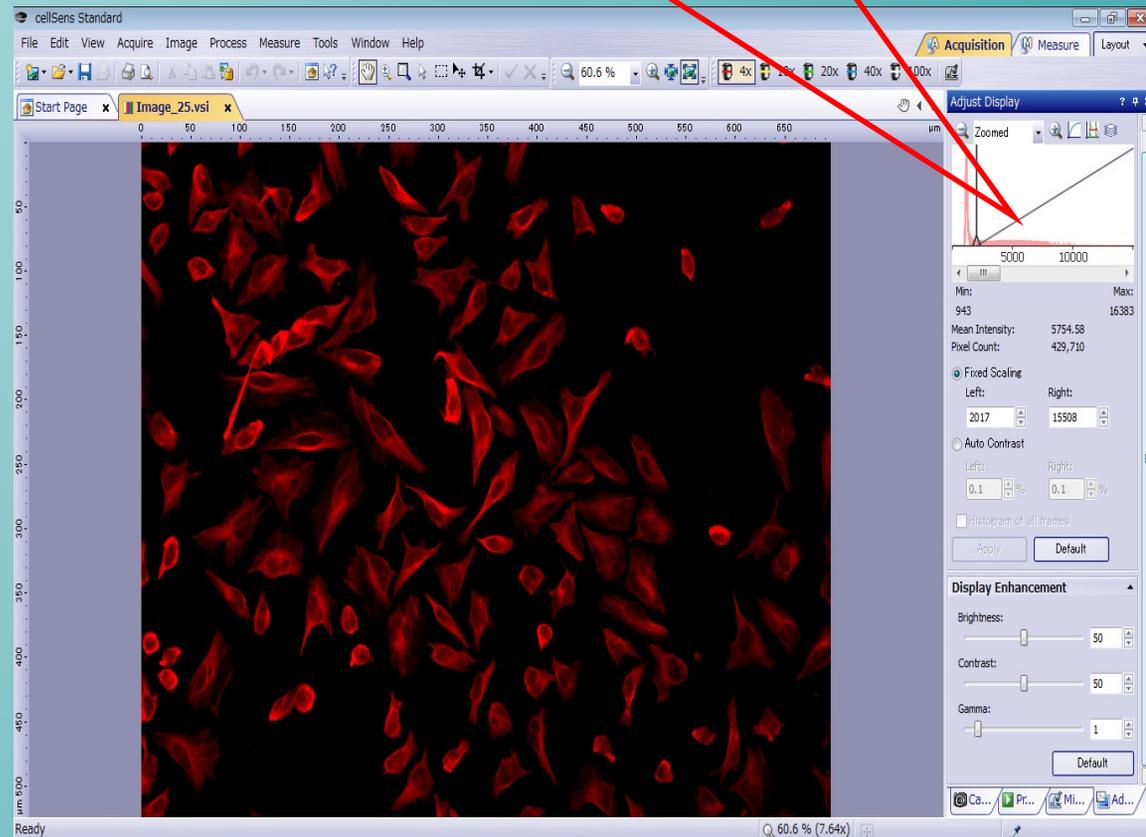
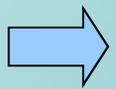
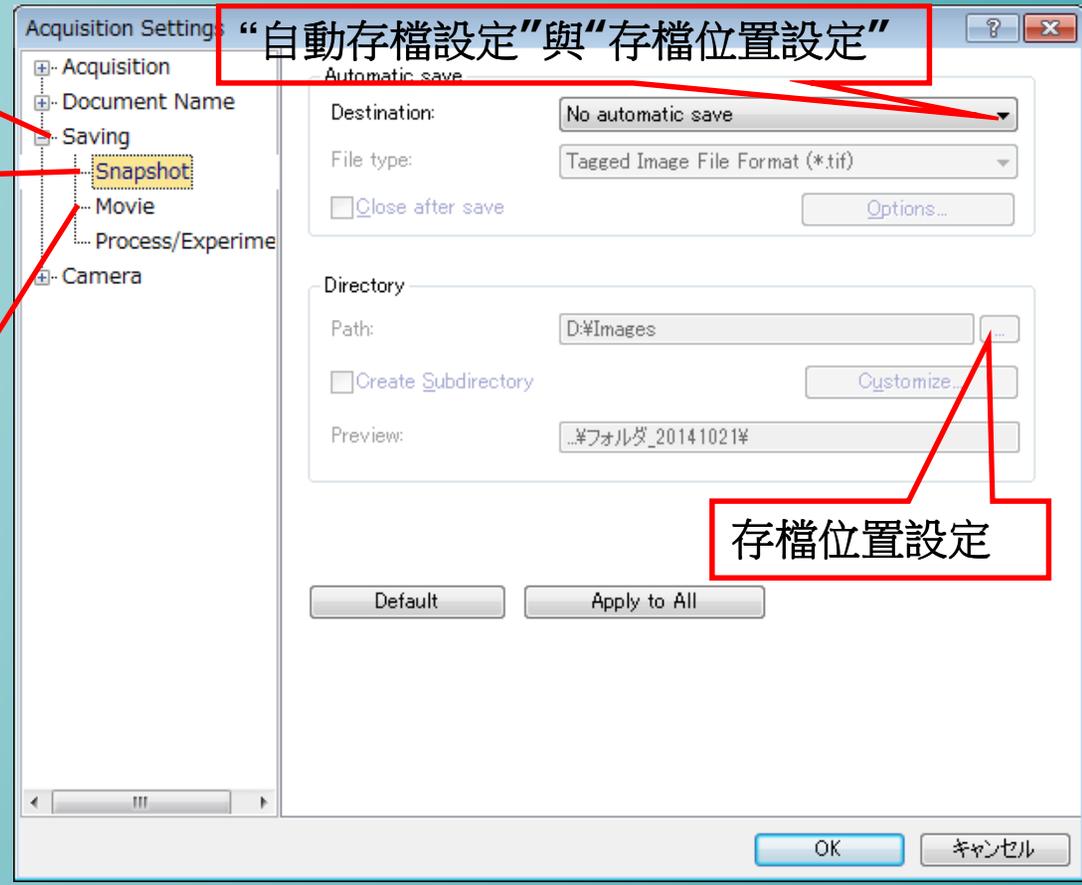
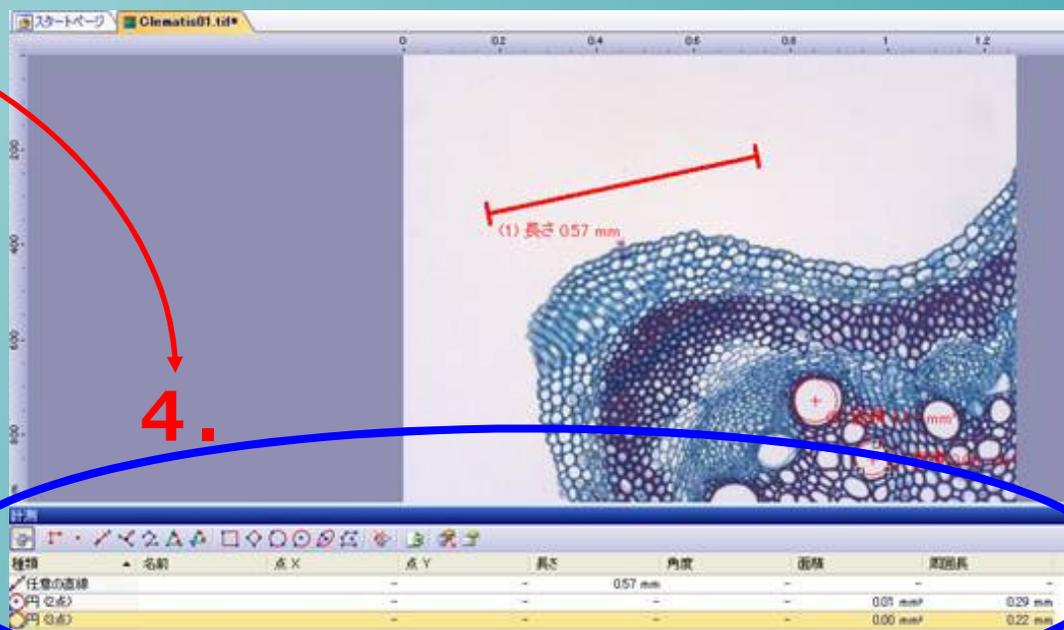
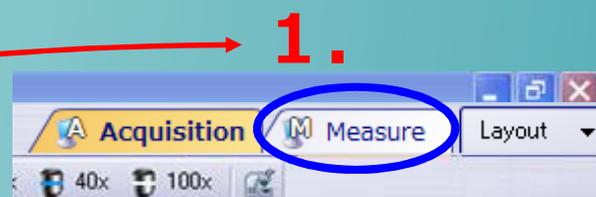
從 **View > Tool Windows > Objectcount** 進入到下面顯示之視窗

The screenshot shows the cellSens Standard interface. The main image displays a field of blue-stained cells with three regions of interest (ROI) labeled 1, 2, and 3. A 'Create Class' dialog box is open, showing a text field for 'Name' and a color selection dropdown. The 'Object Counting' tool window is visible at the bottom, showing a list of classes and a bar chart of results.

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The screenshot displays the CellSens Standard software interface. The 'View' menu is open, and the 'Gallery' option is highlighted. The 'Gallery' tool window is also open, showing a grid of image thumbnails. Red annotations and text provide instructions on how to navigate the interface.

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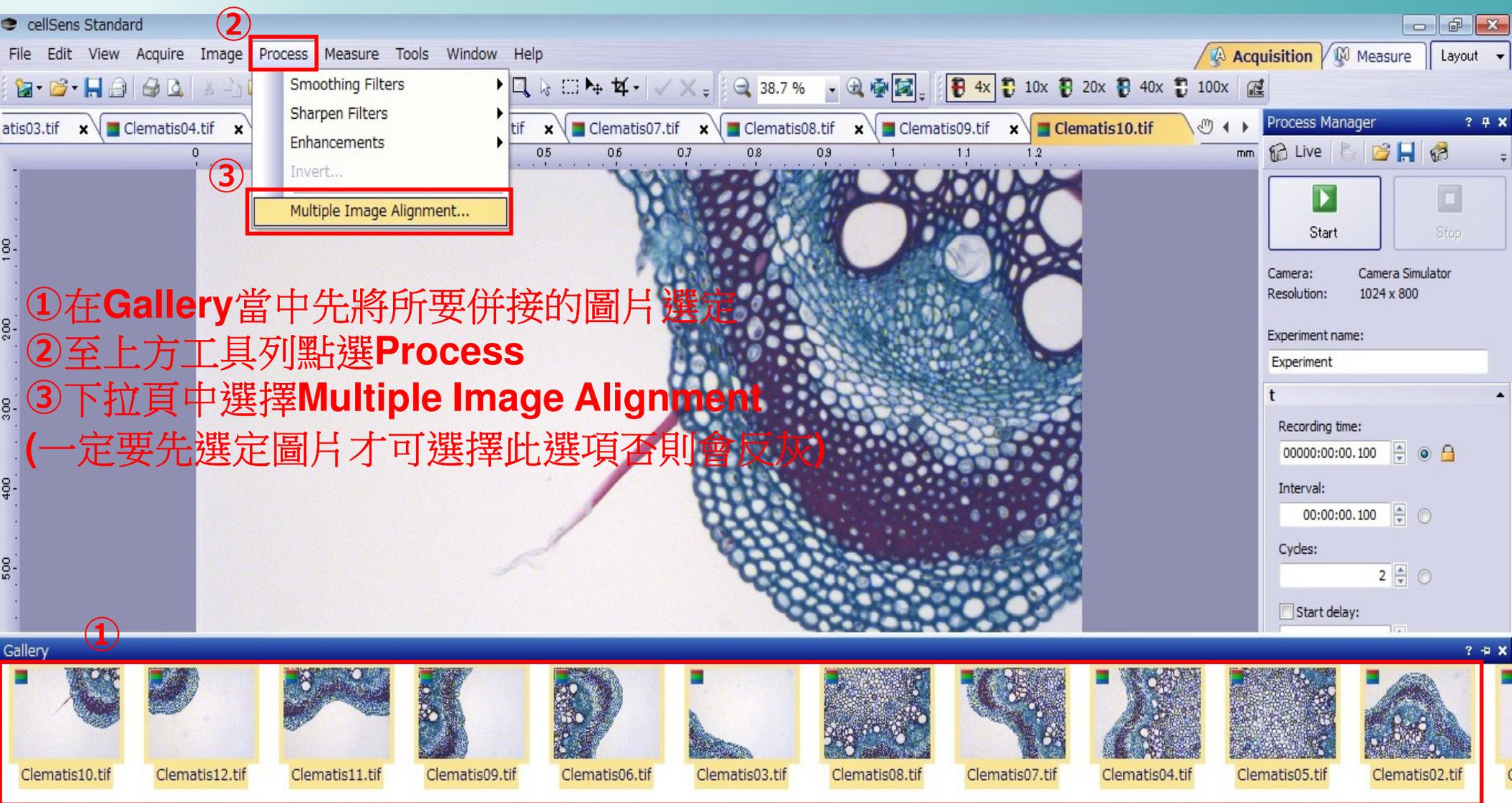
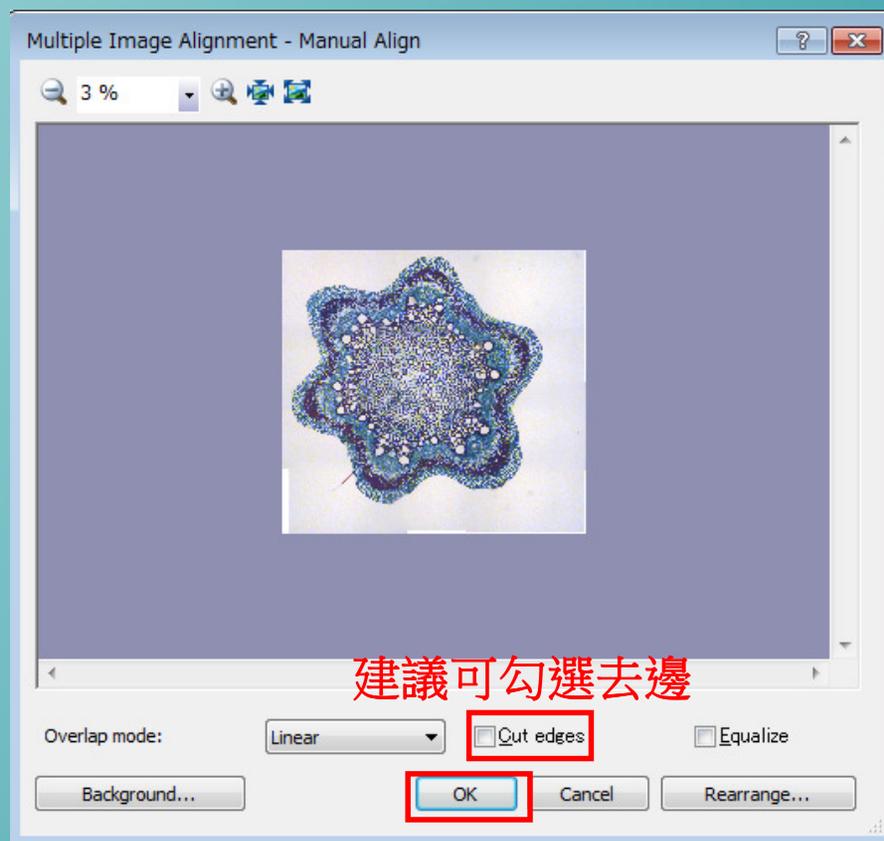
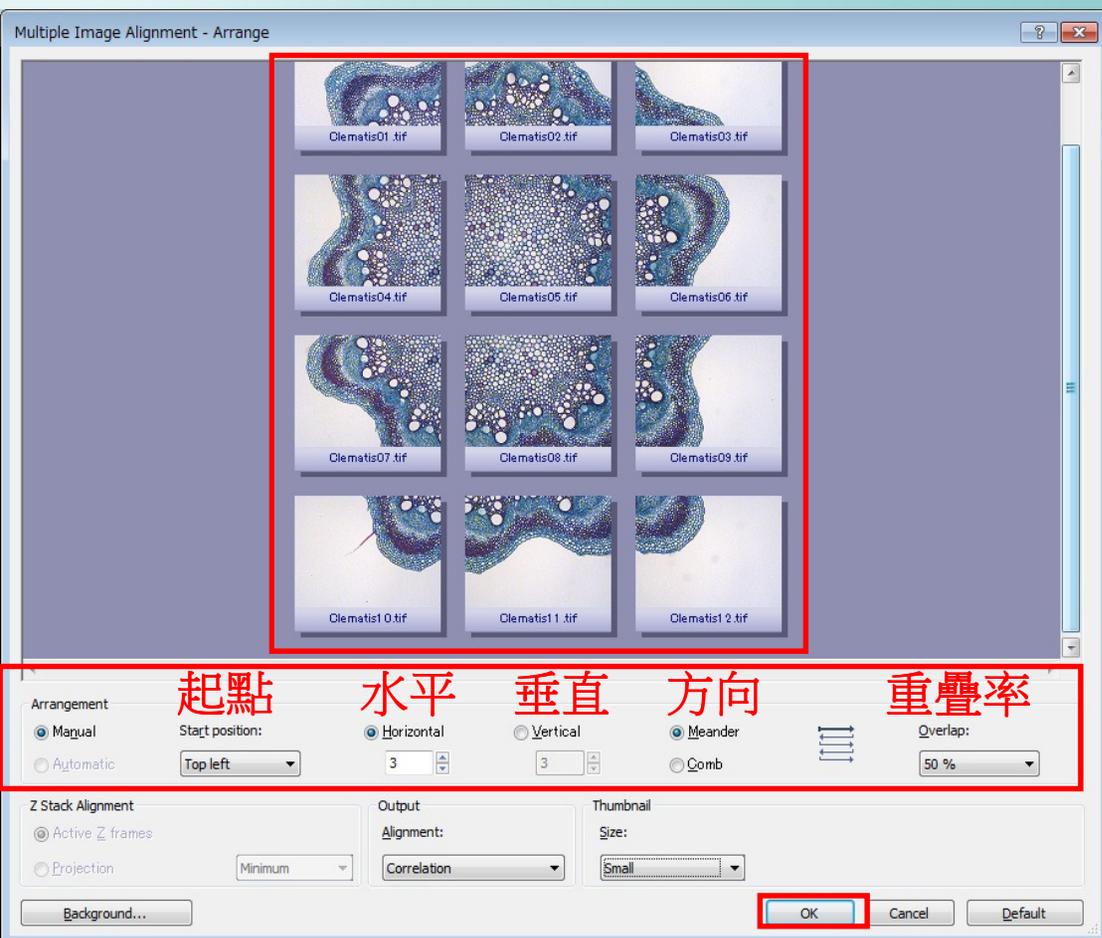
cellSens Standard

File Edit View Acquire Image Process Measure Tools Window

Adjust Display
Pseudo Color...
Mode
Geometry
Burn In Info
Crop...
Calibrate Image...
Change Basic Unit
Separate
Combine Channels...
Combine Color Images...
Duplicate

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cellSens Standard ①

File Edit View Acquire **Image** Process Measure Tools Window Help

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